



MP Biomedicals, LLC

29525 Fountain Parkway
Solon, Ohio 44139

Telephone: 440/337-1200
Toll Free: 800/854-0530
Fax: 440/337-1180
mailto: biotech@mpbio.com
web: <http://www.mpbio.com>

TECHNICAL INFORMATION

Catalog Number: 101076, 193980
Ribonuclease A

Molecular Weight: 13,700

CAS # : 9001-99-4

E.C. # 2.7.7.16

Physical Description: White lyophilized powder

Source: *Bovine pancreas*

Composition: The composition and structure of ribonuclease A has been extensively investigated. The tertiary structure indicates that flexibility in a three-dimensional structure is essential for optimal catalytic activity. ^(2,3,4)

Optimum pH: 7.0 - 7.5 ⁽⁵⁾

Extinction coefficient: $E^{1\%}_{280} = 7.3$

Isoelectric point: pH 9.45 ⁽⁶⁾

Specificity: The enzyme is specific for pyrimidine nucleoside linkages, i.e., those cleaving cyclic 2',3'-pyrimidine nucleotide residues either singly or at the termination of purine nucleotide chains. ⁽⁷⁾

Inhibitors: Ribonuclease is inhibited by heavy metal ions and is competitively inhibited by DNA. The effect of denatured DNA is much greater than that of the native nucleic acid. ⁽⁸⁾

Stability: The enzyme is stable for years stored as a refrigerated dry powder or frozen in solution; refrigerated solutions are stable for weeks. The enzyme aggregates upon lyophilization and storage. ⁽⁹⁾ Special care should be given to handling of the enzyme because of its affinity for glass surfaces ⁽¹⁰⁾

Unit Definition: That amount of enzyme causing the hydrolysis of RNA at a rate such that k (velocity constant) equals unity at 25°C and pH 5.0.

Non-specific protease activity: Not detectable.

Purity: >70%

Solubility: Dissolves readily at 2 mg/ml in analytical grade water to give a clear colorless solution.

Availability:

Catalog Number	Description	Size
101076	Ribonuclease A	100 mg 250 mg 1 g
193980	Ribonuclease A, molecular biology reagent	10 mg 50 mg 250 mg 500 mg 1 g

References:

- Hirs, C., Moore, S., and Stein, W., "Peptides Obtained by tryptic hydrolysis of performic acid-oxidized ribonuclease," *J. Biol. Chem.*, **v. 219**, 623 (1956).
- Kartha, G., Bello, J., and Harker, D., "Tertiary Structure of Ribonuclease," *Nature*, **v. 213**, 862 (1967).
- Lin, M., "The Structural Role of Amino Acid Residues Near the Carboxyl Terminus of Bovine Pancreatic Ribonuclease A," *J. Biol. Chem.*, **v. 245**, 6726 (1970).
- Wang, M., et. al., "Transition-state structure for a conformation change of Ribonuclease," *Bioorg. Chem.*, **v. 4**, 392 (1975).
- Brown, D., and Todd, A., "Cytochrome C peroxidase compound ES is identical with Horseradish Peroxidase Compound I in iron-ligand distances," in *The Nucleic Acids, I*, (Cahrgaff, E., and Davidson, J., eds.), Academic Press, NY, 409 (1955).
- Richards, F., and Wyckoff, H., *The Enzymes, IV, 3rd Ed*, (Boyer, P., ed.), Academic Press, NY, 647 (1971).
- Volkin, E., and Cohn, W., "On the Structure of Ribonucleic Acids. II. The products of Ribonuclease Action," *J. Biol. Chem.*, **v. 205**, 767 (1953).
- Sekine, H., Nakano, E., and Sakaguchi, K., "The interaction of DNA with Pancreatic Ribonuclease A," *Biochim. Biophys. Acta*, **v. 174**, 202 (1969).
- Fruchter, R., and Crestfield, A., "Preparation and Properties of Two Active Forms of Ribonuclease Dimer," *J. Biol. Chem.*, **v. 240**, 3868 (1965).
- Hummel, J., and Anderson, B., "Ribonuclease Adsorption on Glass Surfaces," *Arch. Biochem, Biophys.*, **v. 112**, 443 (1965).